



AS 1.2/E




# ANALYSIS OF INDIGESTION TABLETS (A BACK TITRATION)



### Aim

The aim of this experiment is to find the mass of calcium carbonate in one *Tums* tablet. The manufacturer claims that there is 500 mg in each tablet.

You will do this by a back titration. Back titrations are used for reactions between acids and insoluble bases. The insoluble base is reacted with a known amount of excess acid and then the amount of left over acid is measured by titration with sodium hydroxide solution.

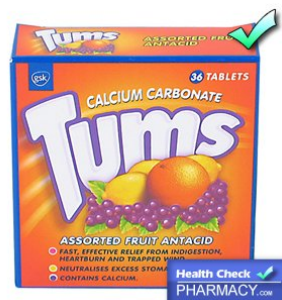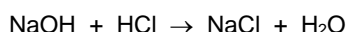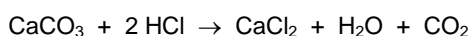

### Safety

	sodium hydroxide hydrochloric acid
--	---------------------------------------

### Method

- 1) Crush up one *Tums* tablet and transfer all of the solid into a volumetric flask.
- 2) Add 25 cm<sup>3</sup> of 0.500 mol dm<sup>-3</sup> hydrochloric acid to the tablet in the volumetric flask. This is an excess.
- 3) Make the mixture up to 250 cm<sup>3</sup> with water in the volumetric flask.
- 4) Titrate 25 cm<sup>3</sup> samples of the stock solution against 0.100 mol dm<sup>-3</sup> sodium hydroxide using phenol red as indicator. Phenol red is yellow in acid and red in alkali. Record your results in a suitable table.

### Analysis

- 5) Calculate the moles of acid added to the indigestion tablet. ....  
.....  
.....
- 6) Calculate the moles of sodium hydroxide that reacted in each titration.  
.....  
.....
- 7) Calculate the total moles of left over acid in the stock solution.  
.....  
.....

8) Calculate the moles of acid that reacted with the calcium carbonate in the tablet.

.....  
.....

9) Calculate the moles of calcium carbonate in the tablet.

.....

10) Calculate the mass of calcium carbonate in the tablet.

.....  
.....

11) Calculate the maximum percentage apparatus error in the final result. Standard errors in apparatus are as follows:

volumetric flask	$\pm 0.1 \text{ cm}^3$
25 cm <sup>3</sup> pipette	$\pm 0.1 \text{ cm}^3$
burette (start & end readings and end point )	$\pm 0.15 \text{ cm}^3$

.....  
.....  
.....  
.....

Evaluation

12) Write down the mean class value for the mass of calcium carbonate in each tablet (we shall assume that the class mean is the correct value).

.....

13) Find the percentage difference between your value and the class mean value.

.....  
.....

14) Comment on the accuracy of your result. ....

.....

15) The manufacturer states that the mass of calcium carbonate in each tablet is 500 mg. Is the manufacturer's figure accurate?

.....  
.....

16) Why is it important that the ingredients in medicines is analysed by independent sources?

.....  
.....

17) If you failed to transfer all of the tablet into the volumetric flask, explain how it would have affected your final value of the mass of calcium carbonate.

.....  
.....